

Quick Guide

Combi 508

Cat. No. : CB-508

Max. RPM : 8,000 rpm
Max. RCF : 8,279 xg

Display Screen

Status of RPM / RCF / Time / Rotor ID / ACC / DEC / Program

Setting Speed


Press once : Setting RPM
Press twice : Setting RCF

Setting Time

Up to < 2 hr
Continuous run (00:00)

View Settings

Verifying the setting parameters during centrifugation

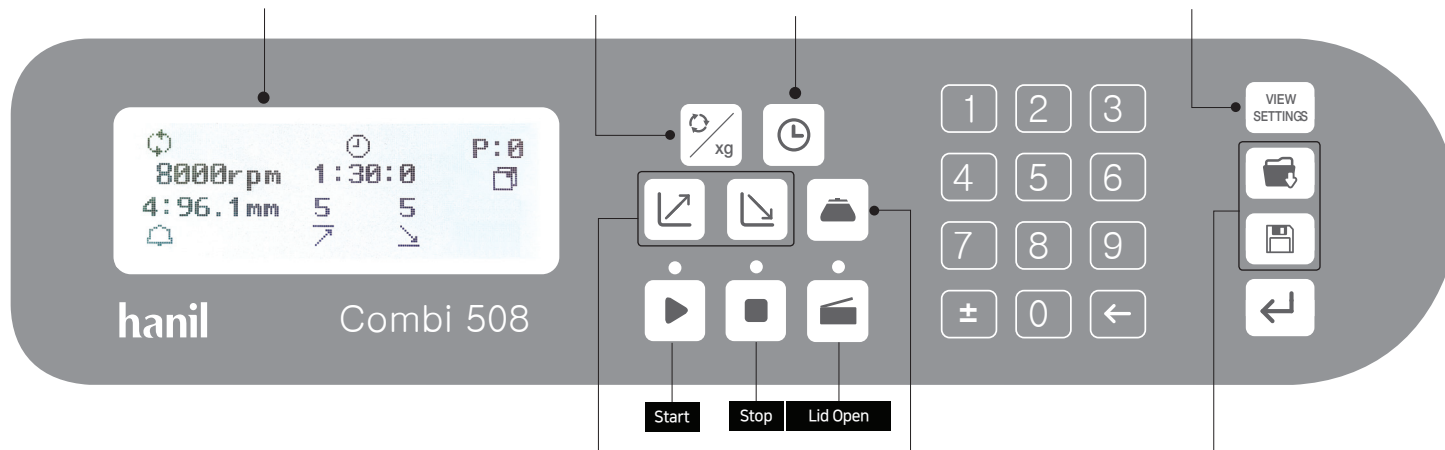

Setting ACC / DEC Rate

- ☒ ACC (0 ~ 9 steps, Max.9)
- ☒ DEC (0 ~ 9 steps, Natural DEC.0)

Setting the Rotor ID manually

When rotor id is not identified automatically, enter the rotor ID. manually

Program Save / Call

- ☒ Saving a program (up to 100)
- ☒ Calling up a program manually

Setting Run Parameters (RPM / RCF / Time / ACC / DEC)



Press a parameter to be changed

Enter the desired value on the keypad

Press the Enter button

Saving / Calling Up a Program




Press the Save or call button

Enter the desired program number on the keypad

Press the Enter button

Cautions for Use

Before Operation

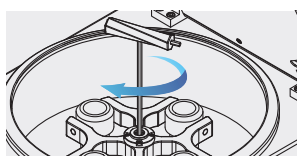

Check the Rotor Locking



Balancing the Centrifuge and Rotor




Load Tubes Symmetrically

After Operation




Keep the Chamber Dry


Maintenance

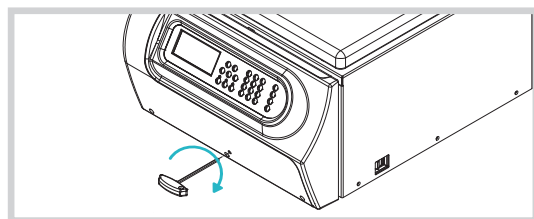

Emergency Lid Open

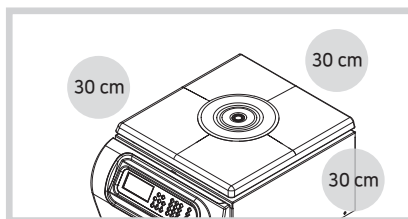

Allow a Safety Clearance of 30 cm

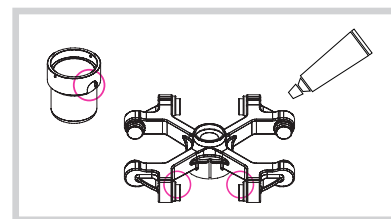

Maintenance of Swing-Out Rotor

hanil

Order & Inquiry | +82-2-3472-0727 | info@ihanol.com | www.ihanol.com
Service | +82-2-3452-8966 | techsupport@ihanol.com




HI-QG-Combi 508(K/E)(Rev.0)

Quick Guide


Combi 508

Cat. No. : CB-508


Max. RPM : 8,000 rpm
Max. RCF : 8,279 xg

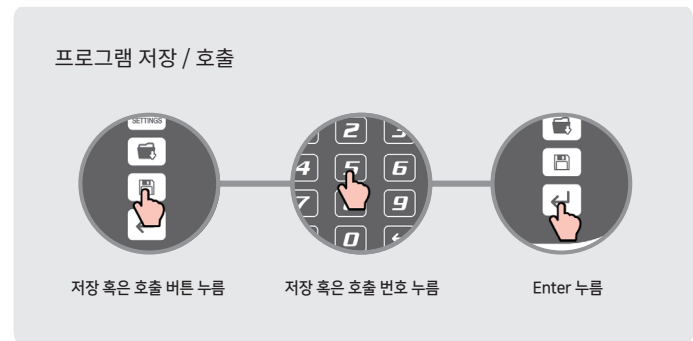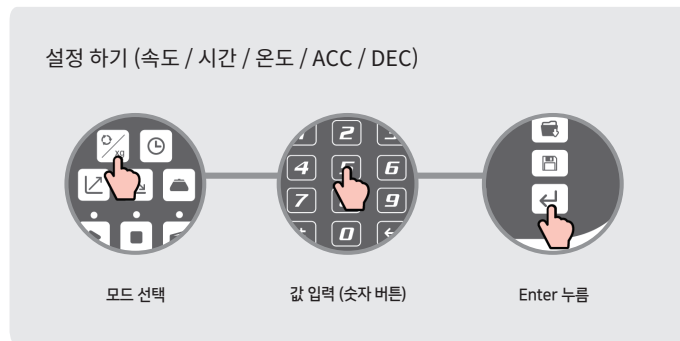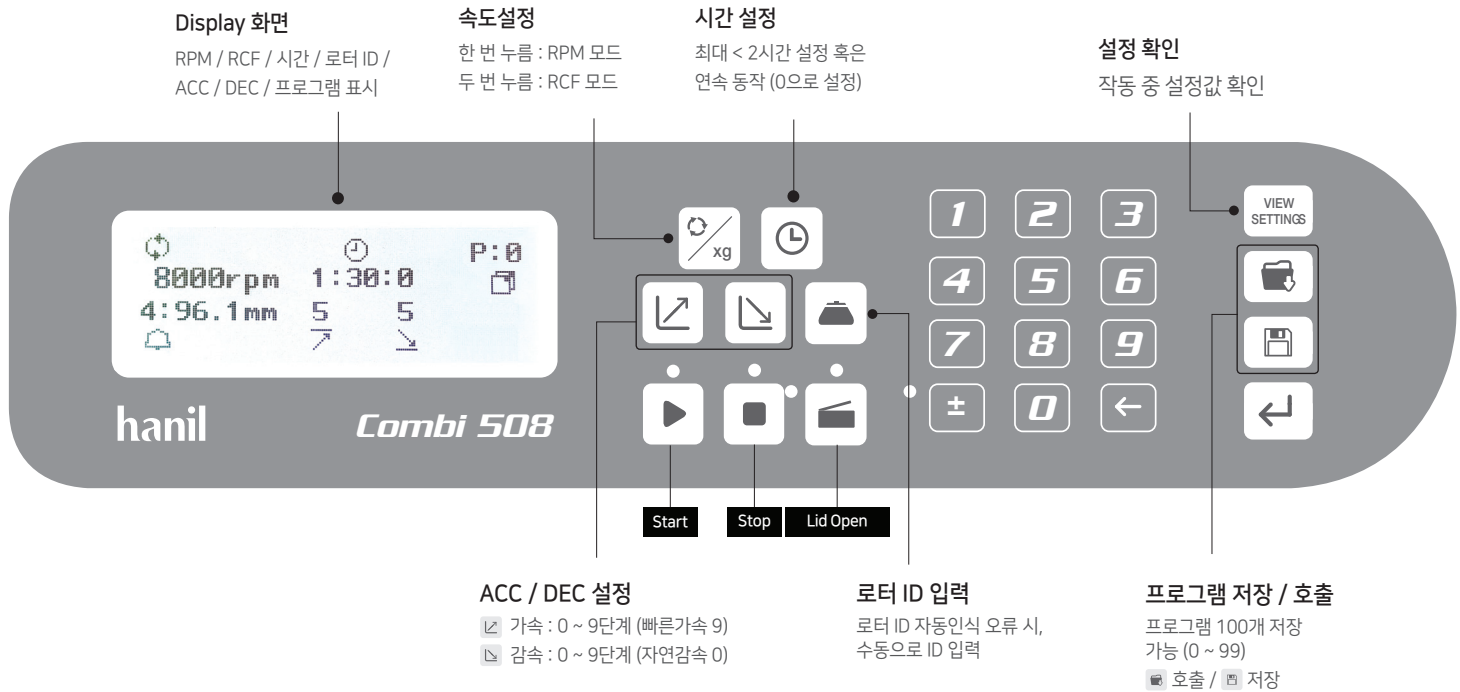


사용 시 주의사항

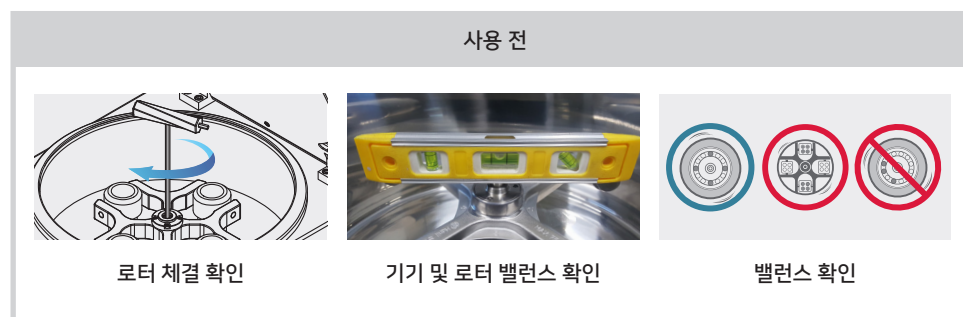


유지관리 주의사항

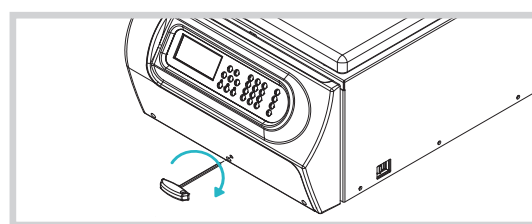

비상 상황 시, 수동 Lid Open

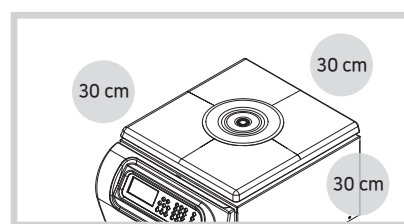

안전 거리 확보

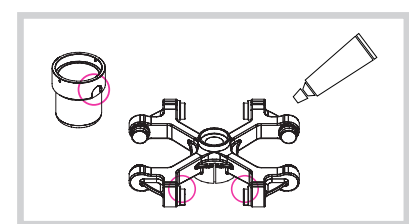

스윙 로터 관리

hanil

제품 문의 | 02-3486-0911 | info@ihanol.com | www.ihanol.com
AS 문의 | 02-3452-8966 | techsupport@ihanol.com